DATASHEET

DETECTORS Photodiode



Photodiode

Vacuum based photodiode



Photek is the world leading manufacturer of vacuum based photodiodes for analysis of ultra-fast optical phenomena in a range of applications.

The vacuum photodiode is a sealed vacuum tube with a much faster response than solid state photodiodes. Vacuum Photodiodes do not contain the MicroChannel Plate electron multipliers used in our Photomultiplier Tubes, resulting in unity gain devices. A benefit of this is the vacuum photodiode's high dynamic range, enabling measurement of extremely high light signals without degradation.

Photek's photodiodes are available in 10 mm, 25 mm and 40 mm formats for use in the UV, visible and NIR spectrum. Photek's photodiodes are the fastest in the world with pulse rise times to 36* ps and pulse FWHM to 55* ps.

Key Attributes

- 10, 25 and 40 mm active areas, other sizes available on request
- > Unity gain
- Wide range of photocathodes including UV, solar blind, visible and NIR response
- > High dynamic range
- > Rise time <36* ps
- Integral 50 ohm output

Applications

- > Analysis of fast optical pulses
- > Cherenkov light detection
- Fluorescence spectroscopy
- > Lidar
- > Particle and nuclear physics
- Measurement of broadband optical phenomena including extreme ultra-violet and vacuum-ultraviolet emission



Options Avaliable


TIME RESPONSE

Time response characteristics are very important when choosing a photodiode (PD) for a particular application, particularly when observing and measuring ultra-fast optical pulses. With the PD010 having the fastest response times, it is the most recommended size. However, 25 mm and 40 mm versions are also available. The pulse rise times and pulse full width half maximum (FWHM) for different sizes of photodiodes are shown below.

Pulse Rise Time				
Diameter (mm)	Minimum (ps)	Typical (ps)	Maximum (ps)	
10	36*	50	60	
25	60	80	100	
40	60	80	100	


Pulse FWHM				
Diameter (mm)	Minimum (ps)	Typical (ps)	Maximum (ps)	
10	55*	60	100	
25	100	150	200	
40	100	150	200	

* Results taken from J. Milnes et al., Recent Developments in Ultra-High Speed and Large Area Photomultiplier Tubes, PoS(ECPD2015)005. The figure below is taken from the paper and shows the influence of the laser pulse width and measurement bandwidth on the observed result:

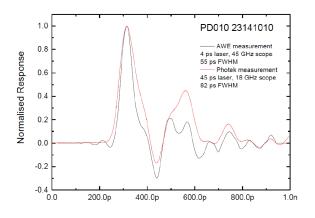


INPUT WINDOW

Photek photodiodes are available with a choice of input window materials. These include MgF₂, fused silica and fibre optic.

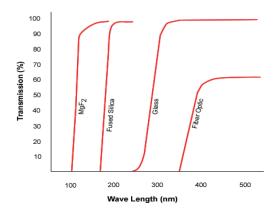

QUANTUM EFFICIENCY CURVES

Photek offer a full range of Gen II photocathodes, these include CsI, Solar Blind, Bialkali, S20 and S25 which demonstrate the broad spectral response that you can achieve as seen below. Please contact the Sales office to discuss your exact requirements.

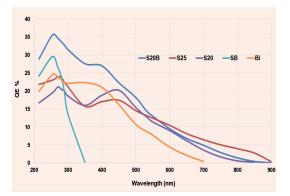

Photek Limited


26 Castleham Road, St Leonards on Sea, East Sussex, TN38 9NS, United Kingdom T: +44 (0)1424 850 555 E: sales@photek.co.uk Registration No. 2641768 England DS041 / rev02 / Aug 21
photek.com




Mechanical

OUTLINE DRAWING


PD010

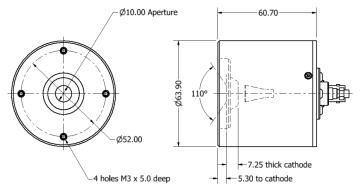

PD025

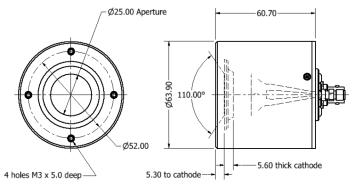

PD040

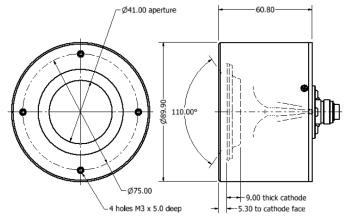

Electrical

-5 kV

Environmental

Operational Limits: -40°C to +45°C Storage: -40°C to +60°C

ENVIRONMENTAL TESTING

For applications where the photodiode is exposed to temperature and shock conditions Photek has the appropriate facilities to offer environmental stress screening. Our vibration system offers shock, sine, random, and sine-on-random testing. Our thermal chamber has a temperature range of -75°C to +160°C and can control humidity from 10% to 98%.

Power Supply

Photek manufactures a range of power supplies for our photodiodes. Our power supplies use the very latest in power supply design and are available in a bench top format.

Pre-Amplifier

Photek manufactures a 2 GHz bandwidth 20 dB gain pre-amplifier, the PA200-10 that is designed to be used with Photek's photodiodes. The PA200-10 is used in close conjunction with the photodiode to amplify the timing pulses for connection by a 50 ohm cable to oscilloscopes or other fast timing electronics.

Photek Limited

26 Castleham Road, St Leonards on Sea, East Sussex, TN38 9NS, United Kingdom T: +44 (0)1424 850 555 E: sales@photek.co.uk Registration No. 2641768 England DS041 / rev02 / Aug 21
photek.com



About Photek

Photek is a specialist manufacturer of vacuum based tubes and camera systems for photon detection.

Our product range includes; Camera Systems, Image Intensifiers, Photomultiplier Tubes, Streak Tubes plus a range of associated electronics.

We are experts in large area and ultra-high speed imaging and advanced photon counting camera systems.

Our continuing success is built upon continuous innovation and product development, and by harnessing and applying knowledge to find solutions for all of our customers' applications.

Photek is accredited to ISO 9001 and ISO 14001.



Contact Us

Our team of specialist engineers and scientists are ready to discuss your application requirements in depth.

- T: +44 (0)1424 850 555
- E: sales@photek.co.uk

Photek Ltd reserves the right to update and improve this document without prior notice.

Photek Limited 26 Castleham Road, St Leonards on Sea, East Sussex, TN38 9NS, United Kingdom T: +44 (0)1424 850 555 E: sales@photek.co.uk Registration No. 2641768 England DS041 / rev02 / Aug 21 photek.com